



# Installing Dry Box Gloves on CBC Glove Extensions with CBC Glove Connecting Kit.

## Step 1

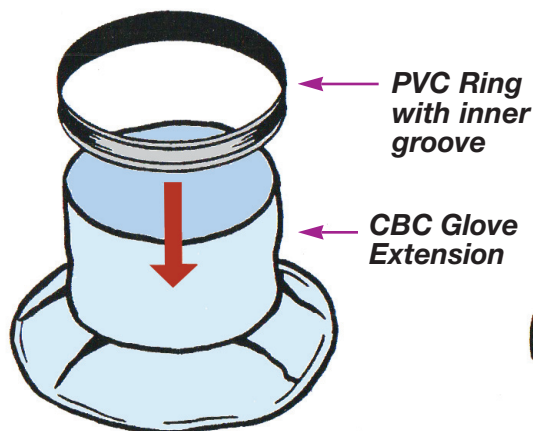

**Step 1.** Place the PVC Ring over the straight portion of the CBC Glove Extension.


## Step 2

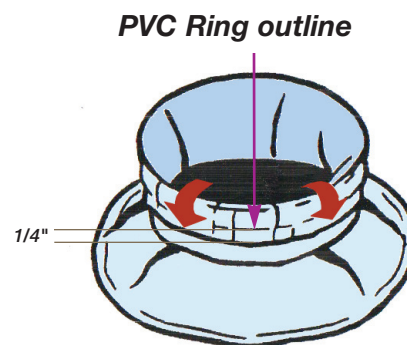

**Step 2.** Fold the straight portion of the CBC Glove Extension over the PVC Ring.

*Note: The edge of the CBC Glove Extension should extend past the PVC Ring by at least 1/4".*


## Step 3

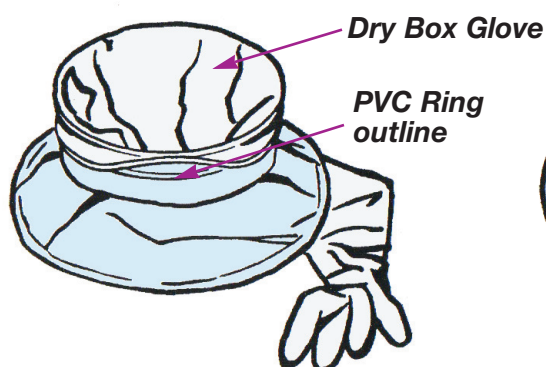

**Step 3.** Insert the Dry Box Glove (with thumb pointing up) through the PVC Glove Extension, leaving enough of the cuff to fold over the PVC Ring outline.

*Note: The bead of the glove should extend beyond the edge to the PVC Ring outline.*


## Step 4

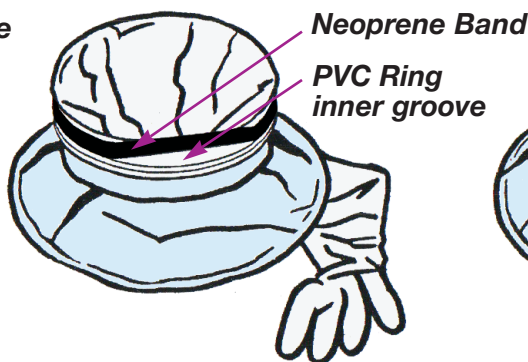

**Step 4.** Place the Neoprene Band over the PVC Ring outline so that it fits within the groove of the PVC ring.

*Note: There will be one thickness of CBC Glove Extension and one thickness of Dry Box Glove between the PVC Ring and the Neoprene Band.*

## Step 5

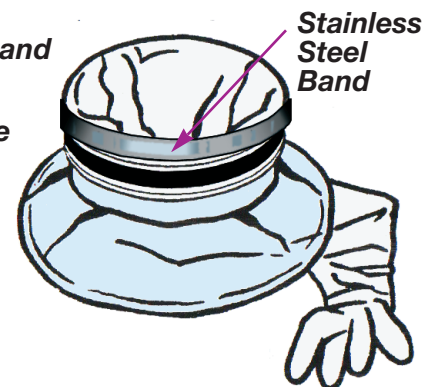

**Step 5.** Place the Stainless Steel Band directly over the Neoprene Band and tighten down.

*Note: The adjusting part of the Stainless Steel Band should be at the top of the CBC Glove Extension and it should be covered with tape or other material to protect the isolator.*

For more information about our CBC Glove Extensions, Dry Box Gloves and other isolator accessories, call **608.273.9661**, or visit our website at: [www.cbcclean.com](http://www.cbcclean.com)



Quality Built to Protect  
Both Research and Researchers

Class Biologically Clean, Ltd • 810 Stewart Street • Madison, WI 53713